

## [CE SNAP SHOT FOR LAUNCH PAD 2.071 +](#)

This program enables you to set up one hand held, copy the settings and then send them to the other hand helds to ensure that all the hand helds are set up the same and properly. It is a good tool and should be used.

First set up one hand held from manufacture settings to working settings, then transfer data to the hand held from the PC. Then set up the system settings in Labrace and then test in manual mode and add the inventory and just be sure everything is working and this is how you want all the hand helds at the location to be set up and working.

### Receive Data from Pocket PC

1. Turn on the hand held and then **Active Sync** it to the office PC.
2. Open **Launch Pad 2.071+**
3. Touch **CE Tools** and then **CE Settings & Database**.
4. Make sure **Active Sync Communication** is checked and hit **Continue**.
5. Touch **Receive Data from Pocket PC**, **wait** and let it do the transfer.
6. Once it is completed then you can check to be sure it worked by doing to the **C drive**, then **program files** and **Launch Pad 2.071+** and then open **CE snapshot** and you should not have two files in this folder "**LabraCE.backup**" and "**LabraCEreg.backup**". These are the files copied in this process.

To set up the other hand helds with this information you do the following steps.

### Send To Pocket PC

1. Turn on the hand held and then **Active Sync** it to the office PC.
2. Open **Launch Pad 2.071+**
3. Touch **CE Tools** and then **CE Settings & Database**.
4. Make sure **Active Sync Communication** is checked and hit **continue**.
5. Touch **Send to Pocket PC**
6. Then you can leave both checked or check the one you want to update to the Hand Held. "**Send Settings**" and or "**Send Database**", hit **OK** and then wait and let it do the transfer.
7. Test to be sure that the data went to the hand held and it is ready for use.

